



BioReady LA Taq

BRPLAT005

- 1 x 500 Units (5 units/ μ l)
- 2 x 10x Reaction Buffer with 15mM MgCl₂ (1ml)
- 1 x 25mM MgCl₂ (1ml)
- 1 x 6x loading dye (500 μ l)

Description

BioReady LA Taq is a thermostable DNA polymerase blend with 3' \rightarrow 5' exonuclease activity.

Storage

-20°C

Definition of Activity

One unit is the amount of enzyme that will incorporate 10 nmol of dNTPs into acid-insoluble product in 30 minutes at 74°C using activated salmon sperm DNA as the template.

Quality Control

This product was purified by HPLC twice.

- | | |
|-------------------------------|-----|
| 1) 5'-3' exonuclease activity | Yes |
| 2) Extra A addition | Yes |
| 3) 3'-5' exonuclease activity | Yes |
| 4) DNase free | Yes |
| 5) RNase free | Yes |
| 6) Protease free | Yes |

Reaction Mixture

10x Reaction Buffer	5 μ l
dNTP Mix (10 mM)	8 μ l
Forward primer (10 μ M)	2 μ l
Reverse primer (10 μ M)	2 μ l
BioReady LA Taq	0.4 μ l
Template DNA	2 ng
ddH ₂ O	to 50 μ l

Thermal Cycler Protocol

94°C	3min	} 35 cycles
94°C	30s	
50-72°C	30s	
72°C	60s/kb	
72°C	5min	

www.bulldog-bio.com | phone: 603-570-4248 | fax: 603-766-0524

One New Hampshire Ave, Suite 125 | Portsmouth, NH 03801

www.bulldog-bio.com | phone: 603-570-4248 | fax: 603-766-0524

One New Hampshire Ave, Suite 125 | Portsmouth, NH 03801