



# Acqua Stain Protein Gel Stain AS001000

1 x 1,000ml bottle

## Description

Acqua Stain is a one-step protein gel dye that requires no pre- or post- gel treatment. Just run your protein gel, add the stain and watch your bands appear.

## Storage

4°C

## Shelf Life

1 year

## Directions

Add enough stain to cover gel completely. Incubate at room temperature according to the times listed below for lowest detectable level in a protein band. Gently agitate if possible.

- |                   |            |
|-------------------|------------|
| 1) 1 ug or more   | 2 minutes  |
| 2) 100 ng to 1 ug | 10 minutes |
| 3) 10ng to 100ng  | 60 minutes |

It is possible to decrease these times by inserting gel and stain into microwave for 20 seconds.

For maximum sensitivity follow staining by a 20 minute gentle agitation in dionized water.

[www.bulldog-bio.com](http://www.bulldog-bio.com) | phone: 603-570-4248 | fax: 603-766-0524

One New Hampshire Ave, Suite 125 | Portsmouth, NH 03801