

# MBP-Trap\_A for Immunoprecipitation of MBP-Fusion Proteins

*Only for research applications, not for diagnostic or therapeutic use.*

**Introduction** Fusion of proteins to *E.coli* Maltose binding protein (MBP) is a common technique to increase solubility and expression level of a protein. Furthermore, MBP- tagged fusion proteins are often used in protein-protein interaction studies. Both approaches require highly specific tools to isolate MBP fusion proteins. MBP-fusion proteins and their interacting factors can be isolated fast and efficiently by immunoprecipitation using the MBP-Trap. MBP-Trap utilizes small recombinant alpaca antibody fragments covalently coupled to the surface of agarose beads.

**Specificity** *E.coli* Maltose binding protein

Content	Reagent	Code	Quantity
	MBP-Trap_A	mbta-20	20 reactions (0.5 ml slurry)
	MBP-Trap_A	mbta-100	100 reactions (2.5 ml slurry)
	MBP-Trap_A	mbta-200	200 reactions (5 ml slurry)
	MBP-Trap_A	mbta-400	400 reactions (10 ml slurry)

**Bead properties** Bead size: ~ 90 µm (cross-linked 4% agarose beads)  
Storage buffer: 20% EtOH  
Binding capacity: 10 µl MBP-Trap\_A slurry binds 0,25 nmol MBP. Binding capacity was determined under optimal conditions with recombinant MBP-fusion protein (74,2 kDa) and will vary depending on the MBP-fusion protein and assay conditions.

**Stability and Storage** Shipped at ambient temperature. Upon receipt store at +4°C.  
Stable for 1 year. Do not freeze.

**Required solutions** **Suggested buffer composition**

Buffer	Composition
Dilution/Wash buffer	10 mM Tris/Cl pH 7.5; 150 mM NaCl; 0.5 mM EDTA
Glycine-elution buffer	200 mM glycine pH 2.5
2 x SDS-sample buffer	120 mM Tris/Cl pH 6.8; 20% glycerol; 4% SDS, 0.04% bromophenol blue; 10% β-mercaptoethanol

Related products	MBP Toolbox	Code
	MBP-binding protein	mbt-250
	MBP-Trap_A Kit	mbtak-20
	Binding control agarose beads	bab-20
	Spin columns	sct-10; sct-20; sct-50

**Support** Please refer to our FAQ section at [www.chromotek.com](http://www.chromotek.com) or contact [support@chromotek.com](mailto:support@chromotek.com)

## Protocol for Immunoprecipitation of MBP-Fusion Proteins using MBP-Trap\_A

We recommend to use 0.5 – 1 mg of protein extract or 1 – 5 µg recombinant protein per Immunoprecipitation.

### Equilibrate beads

1. Vortex MBP-Trap\_A beads and pipette 25 µl bead slurry into 500 µl ice-cold wash/dilution buffer. Centrifuge at 2.500x g for 2 min at +4°C. Discard supernatant and repeat wash twice.

### Bind proteins

2. Add protein extract or recombinant MBP protein diluted in wash/dilution buffer to equilibrated MBP-Trap\_A beads (step 1). If required, save input for immunoblot analysis. Tumble end-over-end for 1 hour at 4°C.

### Wash beads

3. Centrifuge at 2.500x g for 2 min at +4°C. If required, save supernatant for immunoblot analysis. Discard remaining supernatant.
4. Resuspend MBP-Trap\_A beads in 500 µl ice-cold wash/dilution buffer.
5. Centrifuge at 2.500x g for 2 min at +4°C. Discard supernatant and repeat wash twice.  
*optional: Increase salt concentration in the second washing step up to 500 mM.*

### Elute proteins

6. Resuspend MBP-Trap\_A beads in 100 µl 2x SDS-sample buffer.
7. Boil resuspended MBP-Trap\_A beads for 10 min at 95°C to dissociate immunocomplexes from MBP-Trap\_A beads. MBP-Trap\_A beads can be collected by centrifugation at 2.500x g for 2 min at 4°C and SDS-PAGE is performed with the supernatant.
8. *optional instead of steps 8 and 9: elute bound proteins by adding 50 µl 0.2 M glycine pH 2.5 (incubation time: 30 sec under constant mixing) followed by centrifugation. Transfer supernatant to a new tube and add 5 µl 1M Tris base pH 10.4 for neutralization. To increase elution efficiency this step can be repeated.*  
*Note: Efficiency of glycine elution will vary depending on MBP-fusion protein. Glycine elution is not as efficient as SDS sample buffer elution.*

#### Limited Use Label License

The purchase of this product conveys to the buyer the limited, non-transferable right to use the purchased amount of the product and components of the product to perform internal research for the sole benefit of the purchaser. No right to resell this product or any of its components is conveyed expressly or by implication. For information on obtaining additional rights, please contact [licensing@chromotek.com](mailto:licensing@chromotek.com).