



AcquaStain Protein Gel Stain AS001000

1 x 1,000ml bottle

Description

AcquaStain is a one-step protein gel dye that typically requires no pre- or post-treatment. Just run your protein gel, add the stain and watch your bands appear.

Storage

4°C

Shelf Life

1 year

Directions

Add enough stain to cover gel completely. Incubate at room temperature according to the times listed below for lowest detectable level in a protein band. Gently agitate if possible.

- | | |
|-------------------|------------|
| 1) 1 ug or more | 2 minutes |
| 2) 100 ng to 1 ug | 10 minutes |
| 3) 10ng to 100ng | 60 minutes |

It is possible to decrease these times by inserting gel and stain into microwave for 20 seconds.

For maximum sensitivity follow staining by a 20 minute gentle agitation in deionized water.