



FastGene® Agarose Gel Band Cutter

For cutting agarose gel bands

Cat.No.: FG-830, 50 pcs.

STORAGE

Store at room temperature (15-25°C)

FastGene® Agarose Gel Band Cutter is intended for research use only.

2020 March / Rev. 03

CONTENTS

COMPONENTS	2
SPECIFICATIONS	2
PROTOCOL	3
ORDERING INFORMATION	4
CONTACT INFORMATION	4


COMPONENTS

FastGene® Agarose Gel Band Cutter	50 pieces FG-830
FastGene® Gel Band Cutter (size: 6 x 3 mm)	50

SPECIFICATIONS

FastGene® Agarose Gel Band Cutter is a ready-to-use and disposable tool for cutting agarose gel band. Please don't reuse the cutter to avoid cross contamination.

The size of the cutted gel band is 6 mm x 3 mm. The band weight is depending on the percentage and thickness of agarose gel. Please refer to the table below for calculating the weight of gel bands.

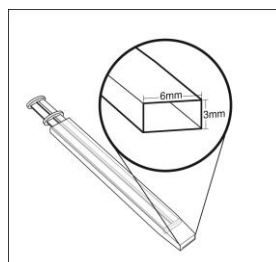


Gel thickness	2% agarose gel	1% agarose gel
1.0 cm	~140 mg	~130 mg
0.8 cm	~120 mg	~110 mg
0.6 cm	~100 mg	~90 mg

PROTOCOL


Agarose Gel Band Cutter

Handling

- Put the gel on the UV box and turn on the power, then use the cutter to cut DNA band directly.
- Remark: Do not press the head of stick when cutting the gel

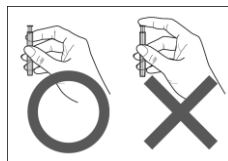

- The cutter can grab DNA bands of the same samples continually and push out gels at one time.

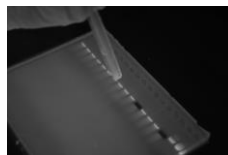

- Push out the gel into microcentrifuge tubes.



ORDERING INFORMATION

Product	Cat.No.
FastGene® Gel/PCR Extraction Kit (100 preps)	FG-91202
FastGene® Gel/PCR Extraction Kit (300 preps)	FG-91302
FastGene® Gel Band Cutter (50)	FG-830
FastGene® Plasmid Mini Kit (100 preps)	FG-90402
FastGene® Plasmid Mini Kit (300 preps)	FG-90502
FastGene® Dye Terminator Removal Kit (50 preps)	FG-9411

CONTACT INFORMATION

EUROPE

Nippon Genetics Europe GmbH
Mariaweilerstraße 28-30
D-52349 Düren
Phone: +49(0)2421-554960
Email: info@nippongenetics.eu
info@nippongenetis.de

For more detailed product information, contact details, questions, or trouble shooting please visit our English website www.nippongenetics.eu or our German website www.nippongenetics.de

JAPAN

Nippon Genetics Co.,Ltd
Head Office
Hongo Tsuna Bldg 3F, 17-9
Hongo 6 Chome, Bunkyo, Tokyo
113-0033 JAPAN
Phone: +81(3)3813-0961
Email: info@genetics-n.co.jp

For more detailed product information, contact details, questions, or trouble shooting please visit Japanese website www.n-genetics.com

FastGene® is a registered trademark of Nippon Genetics Europe GmbH