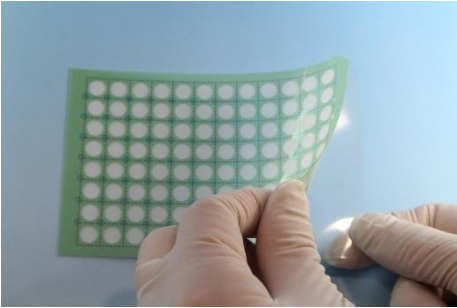
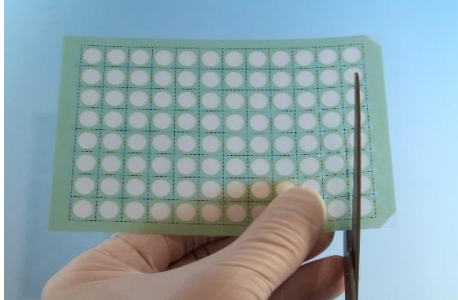
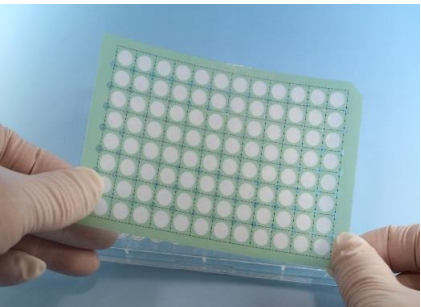
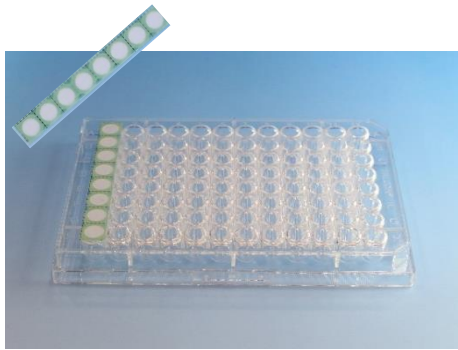
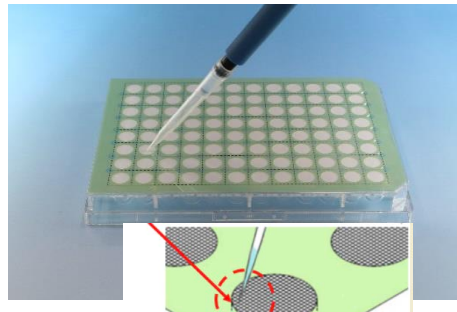
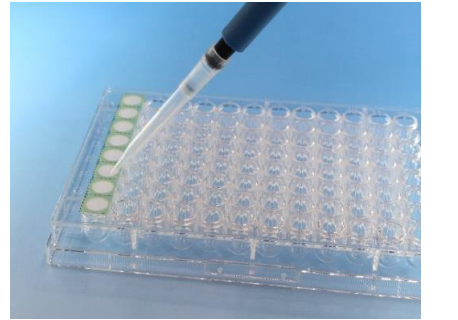


Directions for Bulldog MeshSeal96 Cell Strainers

	96 well use		Custom well number use
①	 <p>Peel off the liner on the back side.</p>	①	 <p>Cut a section off along the perforation line using a scissors.</p>
②	 <p>Place the MeshSeal96 on the microplate centering the filter over the wells. It is possible to remove and replace several times if needed to get the correct fit.</p>	②	 <p>Peel off the liner on the back side and place over desired well(s).</p>
③	 <p>Gently place pipette tip on edge of the mesh/well opening at a 45° angle. Slowly depress the plunger which should carefully push out the cells & media through the filter and along the wall of the side wall of the well.</p>	③	 <p>Gently place pipette tip on edge of the mesh/well opening at a 45° angle. Slowly depress the plunger which should carefully push out the cells & media through the filter and along the wall of the side wall of the well.</p>

*This type of dispensing avoids droplets from forming under the seal which are difficult to then get into the well.